
SVBONY

USB Microscope Digital Camera
User Manual

SV189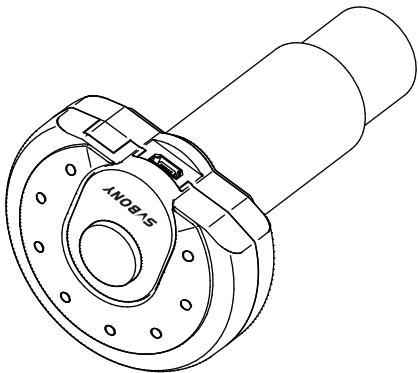

EN

JP

DE

FR

IT

ES

RU

CONTENTS

Product Introduction	1
Installation guidelines	2
Software function introduction	4
Warning	5

◆Product Introduction

Congratulations on your purchase of SVBONY SV189 USB digital microscope camera for microscope. Suitable for microscopes with plug-in standard 23.2/30MM eyepiece interface This manual will briefly introduce you how to use the SV189 USB digital microscope camera.

SVBONY reserves the right to modify this document without notice. Please download the latest version of the user manual pdf file from our official website or print this file.

Please read this manual completely before use. If you have any questions, please feel free to contact us: Facebook: www.facebook.com/svbonny
For software installation instructions and other technical support information, please send us an email: info@svbonny.com

1.Package Including

- ①SV189 digital microscope camera
- ②microscope adapter
- ③1.5m USB 2.0 Cable
- ④CD-ROM
- ⑤User's Manual

2. When using, install the microscope adapter to the camera and tighten it .

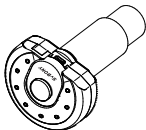
Figure 2. Adapter assembly is complete

3. Specifications

Eyepiece interface size	23.2mm, 30.00mm
Suitable microscope method	plug-in optical microscope eyepiece holder
Photosensitive component	1/4"1.3 million pixel CMOS
Static image resolution	1280*1024, 640*480, 320*240
dynamic image resolution	1280*1024Pixels/24fps
Sensitivity	1.0V/lux-sec@550nm
Pixel size	2.8×2.8um
Spectrum response	400nm---1000nm
Shutter type	ERS (electronic rolling shutter)
Frame rate	16 f/s @ 1280*1024
Image format	JPG/BMP format
White balance	automatic/manual
Interface	USB2.0

◆ Installation guidelines

1. Installation instructions:

PC System Requirements:

Windows(32, 64 bit) XP/7/8/10 /MAC OS10.5 and above / photo booth OS 10.13.3 .

SVBONY official download address:

<https://www.svbony.com/Support/SoftWare-Driver/> Enter the address from the browser to download and install.

Use the attached CD or download "S-EYE_Setup-1.4.4.500" (Figure 3) from the SVBONY official website. After the installation is successful, the shortcut icon "S-EYE" (Figure 4) will be displayed.

Figure 3

S-EYE_Setup-1.4.4.500

Figure 4

S-EYE

2. CD installation:

2.1 Put the CD into the CD-ROM, and the files of the CD will be automatically displayed on the screen.

Note: If the CD-ROM does not run automatically, please double-click "CD-ROM", or right-click "open" under the CD-ROM menu bar

2.2 Double-click the installation file "S-EYE_Setup" to enter the S-EYE Installation Wizard installation wizard. You must accept the user license agreement, so please click: "Accept and Install" to agree to continue the installation . Do not click "Cancel" to exit (Figure 5).

2.3 The default location of the installation directory is "C:\Program Files (x86)\S-EYE", you can click "Browse" to change it to the desired destination folder (Figure 6), check to create a desktop icon , Easy to use and operate.

Figure 5


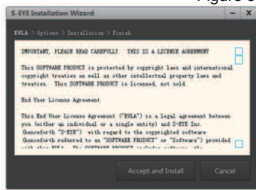
Figure 6

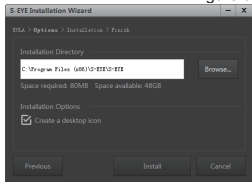2.4 Click "Install" to start the installation and enter the installation process . Do not click "Cancel" to exit (Figure 7). When the S-EYE icon appears on the desktop, the program installation is complete. As shown

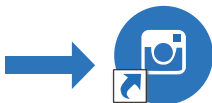
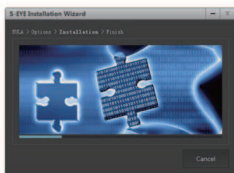
Figure 7

S-EYE

3. Installation/debugging:

3.1 Put the observation specimen into the microscope stage, and fix it with the pressing clip; turn the light source adjustment knob to adjust to the desired light source position, turn the low-magnification objective lens to the positioning point, and move the horizontal and vertical adjustment on the stage Ruler, make the observation specimen to the center position; fine-tune the focus adjustment wheel to obtain a clear imaging effect, take out the eyepiece and extender on the microscope, and insert the SV189 microscope digital camera into the eyepiece tube (Figure 8);

3.2 Connect the SV189 microscope digital camera to the computer. Note: Before connecting the USB camera to the computer, please make sure that the USB2.0 interface on the computer can work normally.

Connect the USB2.0 cable of SV189 to the computer, Open "S-EYE", click the camera name "1.3M Camera" in the camera list in the upper right corner of the software interface, the blue indicator light of the SV189 microscope digital camera will light up and enter the power-on state; SV189 imaging is displayed on the screen (Figure 9) The imaging display

Figure 8

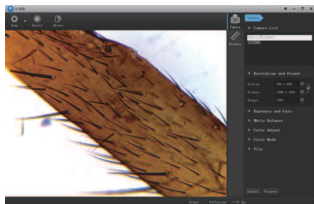
Figure 9

effect can be achieved by manually fine-tuning the focus adjustment wheel to achieve a clearer effect.

◆ Software function introduction

1. Software interface operation: After entering the "S-EYE" software interface, In the "Camera" mode, Snap /Record/Browse functions can be operated, and Resolution and Format/Exposure and Gain/White Balance/Color Adjust/Color Mode/Flip can be set individually (Figure 10)

2. In the "Measure" mode, Use the measuring tool to measure specimen Width/ Height /Radius/ Area /Length /Angle/Distance and other information (Figure 10)

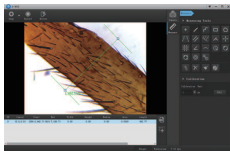
Figure 10

3. Click the "Output" function to view the measurement data list and export the measurement record data to a file.

Click the " " Capture preview image to a file, including measurement tools on the image.


Click the " " Export measurement data to csv file.

◆ Introduction to measurement toolse

	Point		Line
	Broken line		Rectangle
	Polygon		Distance between two lines
	Parallel line		Distance to the first line
	Perpendicular		Crosshair
	Two Crosshairs		Angle
	Arc		Circle
	Two points circle		Three points circle
	Concentric circles		The distance between the two circles
	Move		Remove
	Remove all measurement tools		Set the color of the measurement tool

4. Click the "Edit" function to calibrate multiple measurement data of the specimen, and "Backup" can backup and save the calibration data as a file (Figure 11).

Figure 11

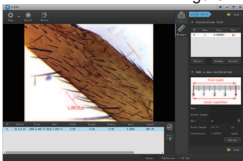◆ Warning

Before using this device, read this guide which contains important operating instructions for safe usage and control for compliance with applicable standards and regulations.

FCC Requirements:

•Products authorized under Part 15 using SDoC or Certification require a label containing one of the following compliance statements

(1)Receivers associated with licensed device service operations:

This device complies with part 15 of the FCC Rules. Operation is subject to the condition that this device does not cause harmful interference.

(2)Stand-alone cable input selector switch:

This device complies with part 15 of the FCC Rules for use with cable television service.

(3)All other devices:

•This device complies with part 15 of the FCC Rules. Operation is subject to the following two conditions:

(1) This device may not cause harmful interference, and

(2) this device must accept any interference received, including interference that may cause undesired operation.

CE Requirements:

•(Simple EU declaration of conformity) Hong Kong Svbon Technology Co.,Ltd declares that the equipment type is in compliance with the essential requirements and other relevant provisions of RED Directive 2014/30/EU and the ROHS Directive 2011/65/EU and the WEEE Directive 2012/19/EU; the full text of the EU declaration of conformity is available at the following internet address: www.svbony.com.

•Disposal

The crossed-out wheeled-bin symbol on your product, literature, or packaging reminds you that in the European Union, all electrical and electronic products, batteries, and accumulators (rechargeable batteries) must be taken to designated collection locations at the end of their working life. Do not dispose of these products as unsorted municipal waste. Dispose of them according to the laws in your area.

IC Requirements:

CAN ICES-3(B)/NMB-3(B)

Avoid Choking Hazard

Small Parts. Not for children under 3 years.

Approved Accessories

WARNING

- This device meets the regulatory standards when used with the Svbonny accessories supplied or designated for the product.
- For a list of Svbonny-approved accessories for your item, visit the following website: <http://www.Svbonny.com>

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EU Importer:

Name: Germany Retevis Technology GmbH

Address: Uetzenacker 29,38176 wendeburg

Warranty Card

Product Model	
Purchasing Date	
Defect Reason	
Dealer Name	
Telephone	
User's Name	
User's Address	
User's Email	

Remarks:

1. This guarantee card should be kept by the user, no replacement if lost.
2. Most new products carry a one-year manufacturer's warranty from the date of purchase.
3. The user can get warranty and after-sales service as below:
 - Contact the seller where you buy.
4. For warranty service, you will need to provide a receipt proof of purchase from the actual seller for verification

Exclusions from Warranty Coverage:

1. To any product damaged by accident.
2. In the event of misuse or abuse of the product or as a result of unauthorized alterations or repairs.
3. If the serial number has been altered, defaced, or removed.

MADE IN CHINA

Hong Kong Svbonny Technology Co.,Ltd

Add: Unit B, 5th Floor, Gallo Commercial Building,
114-118 Lockhart Road, Wanchai, Hong Kong

Facebook: [facebook.com/svbony](https://www.facebook.com/svbony)

E-mail: info@svbony.com

Web: www.svbony.com

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