
SVBONY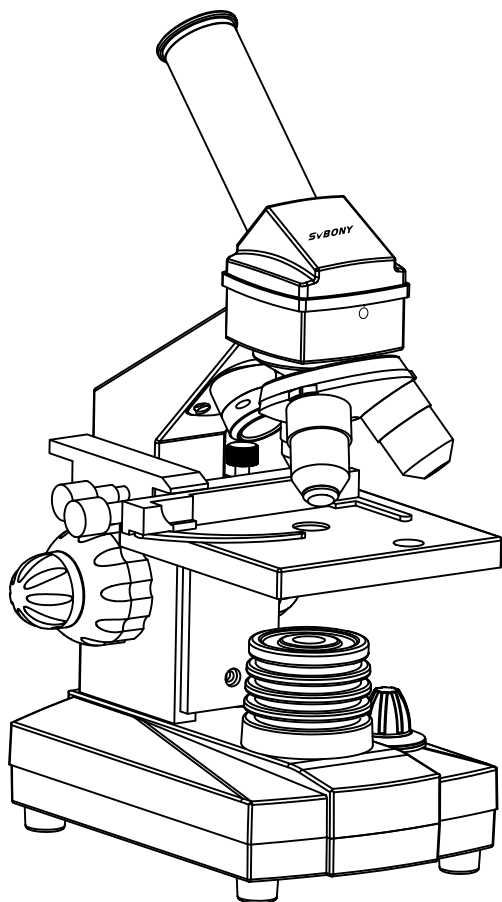

EN

JP

DE

FR

IT

ES

SV601

PREFACE

Thank you very much for your purchase of a SVBONY microscope
This manual applies to SV601. Read this manual before using the SV601 microscope.

Setting Up

1. Remove the Styrofoam container from the carton.
 2. Carefully remove the microscope and other parts from the container and set them on a table, desk, or other flat surface.
 3. Remove eyepieces from the plastic bags.
 4. Insert the 10x eyepiece into the eyepiece tube
- You are now ready to use your microscope for looking at specimen


CAUTION

Please set the microscope on stable table, desk, or other flat surface. Away from edge.

●CARE & MAINTENANCE

1. Never leave the SV601 microscope close to fire, or hot places
2. When you are done using your microscope, remove any specimens left on the stage
3. Turn off the top and bottom illuminators when you are done using the microscope.
4. If you will not use the microscope for a long time (30 days or more), remove the batteries.
5. Always place the dust cap over the eyepiece when not in use or when being stored.
6. Store the microscope in a clean, dry place.
7. Clean the outside surfaces with a damp cloth.
8. Never clean optical surfaces with cloth or paper towels as they can scratch optical surfaces easily.
9. To remove fingerprints from optical surfaces, use a lens cleaning agent and lens tissue available at most photo outlets.
10. Never disassemble or clean internal optical surfaces. This should only be done by qualified technicians at the factory or other authorized repair facilities.

OVERVIEW

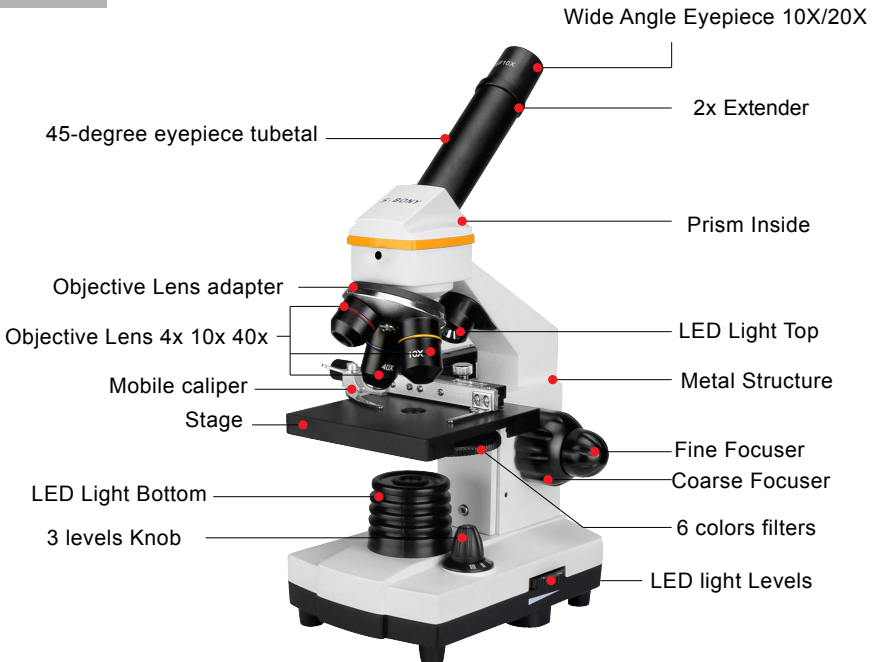List and View of Accessory

SPECIFICATIONS

Model	SV601
Optical System	Finite Biological Microscope
Eyepiece	WF10X WF20X
Objective	4X 10X S40X
Extender	2.0X
Working Stage	90mmX90mm Square
	Moving Scale
	Disc Filters
Adjustment	Coarse Focuser & Fine Focuser (dual knobs)
Illuminator — Top	LED Lights
Illuminator — Bottom	LED Lights , Uses 2AA batteries (user supplied)
Accessories	Full Set (Accessory Chart)
Packing	Box
Dimension	155x110x290mm
Weight	1.5kg

SETTING UP & OPERATION

- ① Remove the microscope, and insert WF10X eyepiece
- ② Insert the battery or connect to electricity, turn on the lights. Make the light in view
- ③ Put the slides on the working stage, fixed by moving scale, adjust object position to the middle of light

- ④ Adjust the objectives to the low position, then move it up slowly by focuser. Adjust fine focuser can get better view

⑤ Move the objectives adaptor, change the magnifications from Min. to Max.

⑥ The filters under working stage can make the vision in different colors.

Precautions

1. Please carefully place the sample in the center of the platform when you use it.
2. Rotate the knurled ring on the objective lens housing to the 4x position. It is best to begin viewing a specimen at low power, then switch to higher power as needed.
3. Look through the eyepieces and gradually turn the focus knob until the specimen comes into focus.
4. If there is not enough travel in the focuser mechanism to reach focus, it may be necessary to lower the stage plate to the down side. and make it travel full distance slowly, until the image becomes sharp.
5. If higher power is desired, rotate the objective lens housing to the 3x position and refocus until the specimen becomes sharp. You can also replace the WF 10x eyepieces with the WF 20x eyepieces to obtain two additional magnifications powers.

- Do not view the sun or other intense light with your products or even with the naked eye; Otherwise there may be permanent eye damage.
- Kids need to be used under the guardianship of adults to avoid the unnecessary damage.
- Do not put the plastic package a place your kids can reach. If not, the kids may put the plastic into the mouth and suffer from suffocating.

Care and Clean:

The item do not need routine maintenance other than making sure that the objective lenses and eyepieces are kept clean. Dirty objectives and/or eyepieces mean less light transmission and loss of brightness as well as unsharp images. Keep your optics clean! When not using your item, store those in the case provided.

Usage Notice:

1. Store the product in a ventilated, dry, clean place or put it in a dry box.
2. If there are dust and other dirt on the optical lens, use a professional cleaning kit to clean it. Do not wipe it directly with a cloth and a hard brush.
3. Please do not disassemble or repair the product yourself. If you need to repair it, please contact our professional after-sales department.
4. Do not use non-waterproof products in wet environments, use waterproof products in wet environments, and dry them after use.
5. When viewing indoors, please do not close the window and watch it through the window, which will reduce the product observation effect.

Avoid Choking Hazard

Small Parts. Not for children under 10 years.

Approved Accessories

- This item meets the regulatory standards when used with the Svbonny accessories supplied or designated for the product.
- For a list of Svbonny-approved accessories for your item, visit the following website: <http://www.Svbonny.com>

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User's Email: _____

Remarks:

1. This guarantee card should be kept by the user, no replacement if lost.
2. Most new products carry a one-year manufacturer's warranty from the date of purchase.
3. The user can get warranty and after-sales service as below:
 - Contact the seller where you buy.
4. For warranty service, you will need to provide a receipt proof of purchase from the actual seller for verification


Exclusions from Warranty Coverage:

1. To any product damaged by accident.
2. In the event of misuse or abuse of the product or as a result of unauthorized alterations or repairs.
3. If the serial number has been altered, defaced, or removed.

EU Importer:

Name: Germany Retevis Technology GmbH

Address: Uetzenacker 29,38176 wendeburg

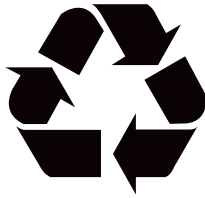MADE IN CHINA

Hong Kong Svbonny Technology Co.,Ltd

Add: Unit B, 5th Floor, Gallo Commercial Building, 114-118
Lockhart Road, Wanchai, Hong Kong

Facebook: [facebook.com/svbonny](https://www.facebook.com/svbonny)

E-mail: info@svbonny.com

Web: www.svbonny.com

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