
SVBONY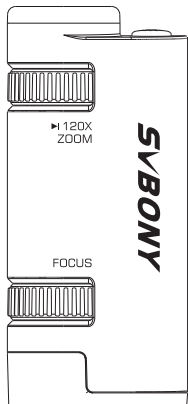

Hong Kong Svbon Technology Co.,Ltd

Add: Unit B, 5th Floor, Gallo Commercial Building, 114-118
Lockhart Road, Wanchai, Hong Kong

Facebook: facebook.com/svbon

E-mail: info@svbon.com

Web: www.svbon.com

User Manual
SV603 Portable Microscope

EN DE FR IT ES

Contents

EN (P01 to P09)

Description	01
Microscope Accessories	02
Operation	02
Specifications	04
Warning	04
Product Protection and Maintenance .	05
Battery Warning	05
Disclaimer	07
Warning	08

DE (P10 to P19)

FR (P20 to P29)

IT (P30 to P38)

ES (P39 to P48)



Caution

Read this user manual before using the SV603 Portable microscope.

Description

Thanks very much for buying SVBONY SV603 portable microscope (SV603). Before using the SV603, please read this user manual patiently, to know more about the functions and operation skills. So that you would have more pleasant experience.

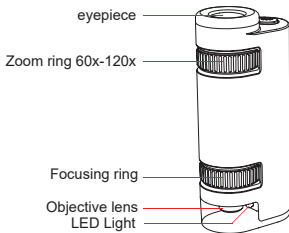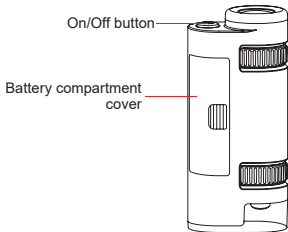Microscope Accessories

1x bag

2x Specimen

(Bee leg slices & Onionskin slices)

1x AAA battery

Operation

Install battery

Due to the SV603 portable microscope has high magnification power (60x-120x), if the microscope illumination is not turned on, it may be low light and affect the observation effect, so it is recommended that you install the battery and turn on the LED lighting.

1. Remove the battery compartment cover.
2. Insert 1 AAA battery.
3. Install the battery compartment cover back to its original position.
4. Turn on the LED light.


Note: The LED light has 3 levels: low light, strong light, and off.

Users can choose different light intensities according to their needs.

Operation

The SV603 portable microscope has 2 Specimens in the package. Choose one specimen beginning the observation. Do as followings:

1. Place the specimen facing up and horizontally on a flat table. Put a piece of white paper with a smooth surface under the slice for a better experience.
 2. Aim the microscope objective lens at the specimen and gently place it on the section to turn on the LED light.
 3. Align one eye to the eyepiece for observation. If the specimen is not observed, you can slowly follow the deviation direction. Slowly adjust to the proper position.
 4. Adjust the focus ring, It is recommended to start observation from the lowest magnification 60x. Turn the focus ring until the image is clear.
- Note: Do not forcefully squeeze the specimen section during observation to avoid damage to the product or section.
5. Please turn off the LED lights in time after the observation.

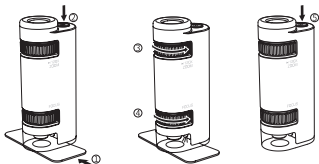Specifications

Product Name: SVBONY SV603 Portable
Microscope

Eye Relief: 10~14mm

Model Number: SV603

Liner Field of View: 1.3~1.9mm

Magnification Power: 60x-120x

Illumination: LED light

Battery Type: AAA 1.5V Battery

Lighting switch: On/Off Button

Length/width/height: 40x25x89mm

Net Weight: 33.5g

Warning

1. Never use SV603 to directly watch the sun or other strong luminous objects, to avoid causing damage to your eyes into permanent damage.
2. Never look directly at the built-in LED luminous body of this product at close range to avoid discomfort to your eyes or pain.
3. When not using the product for a long time, please remove the battery and keep it in a safe place to avoid battery leakage and damage the product.
4. Using this product for a long time will cause the product to heat up, which is normal.
5. This product contains tiny parts. For your child's safety, when provided to your child, parents must be accompanied during the whole process. After use, please keep the product out of children's reach.

Product Protection and Maintenance

If your SV603 microscope is used and kept properly, it can provide you with years of reliable service.

1. Please protect the microscope, keep it away from impact, use it within the limits of moving parts, and do not use brute force.
2. Please protect the microscope lens, and put it into the protective bag when not in use to avoid collision.
3. Please try to keep your microscope in a dry and cool place.
4. When storing for a long time, please put the microscope in a plastic bag or sealed container with desiccant inside.
5. Please do not put the microscope in a car under the hot sun or close to any heat source, these can cause damage.
6. Please clean off dust, dirt or moisture as soon as possible after use, so as not to enter the microscope or enter the mobile phone. Moving parts, causing unpredictable damage.

Battery warning

Please read the instructions and warnings completely before using this equipment. Misuse of this product may cause product damage, overheating, toxic gas, fire or explosion. If harm is caused to you (the buyer), Svbonny (including the manufacturer) is exempt.

- Never store the battery in high temperature environments, including strong sunlight. Never put the battery in the fire or other overheated environment.

- It is forbidden to drop, bump, wear or other actions that damage the battery. If the battery happens any of the following damages, such as dents, bulges, perforations, tearing, deformation or corrosion, etc., no matter how it is caused, please stop using it immediately and contact the manufacturer or hand it to your local battery recycling center for disposal.
- Never disassemble the battery or repair or modify the battery in any way.
- Do not get the battery wet or immerse it in liquid. Keep the battery dry at all times.
- Do not replace any part of the battery (replace the entire battery).
- The battery is designed for adult use. If the purchaser intends to use it for minors, the purchaser agrees to provide detailed instructions and warnings before use. If this is not done, the purchaser shall bear full responsibility because he agrees to the manufacturer's exemption clause for unintended use/abuse by minors.
- All batteries have undergone strict quality inspection. If you find that the battery overheats, smells, if there is abnormal shape, abrasion, fracture, or abnormal phenomena, stop using all batteries immediately and contact the manufacturer.
- The standard battery is a non-rechargeable battery and must not be charged. If not used for a long time, please take it out and store separately.
- Do not put the battery in the trash. Under the terms of environmental laws and regulations, it is illegal to discard used batteries directly in

ordinary trash cans. Please take the used batteries to the local battery recycling center for disposal.

Disclaimer

This battery should only be used with appropriate equipment. Please refer to the packaging of the device to determine if compatible with your specific device.

The manufacturer does not assume any responsibility for any damage to any equipment caused by the use of this product.

The manufacturer does not assume any responsibility in any form, including any damages to you or any third party that you or any third party may cause due to use, whether intentional or unintentional, or misuse of this battery in combination with any equipment or accessories.

The manufacturer will not be responsible for any losses that may be caused by you or any third party misusing this battery as described above.

WARNING

- Do not view the sun or other intense light with your products or even with the naked eye; Otherwise there may be permanent eye damage.
- Kids need to be used under the guardianship of adults to avoid the unnecessary damage.
- Do not put the plastic package a place your kids can reach. If not, the kids may put the plastic into the mouth and suffer from suffocating.

Care and Clean:

The item do not need routine maintenance other than making sure that the objective lenses and eyepieces are kept clean. Dirty objectives and/or eyepieces mean less light transmission and loss of brightness as well as unsharp images. Keep your optics clean! When not using your item, store those in the case provided.

Usage Notice:

1. Store the product in a ventilated, dry, clean place or put it in a dry box.
2. If there are dust and other dirt on the optical lens, use a professional cleaning kit to clean it. Do not wipe it directly with a cloth and a hard brush.
3. Please do not disassemble or repair the product yourself. If you need to repair it, please contact our professional after-sales department.
4. Do not use non-waterproof products in wet environments, use waterproof products in wet environments, and dry them after use.

5. When viewing indoors, please do not close the window and watch it through the window, which will reduce the product observation effect.

Avoid Choking Hazard

Small Parts. Not for children under 3 years.

Approved Accessories

WARNING

- This item meets the regulatory standards when used with the Svbonny accessories supplied or designated for the product.
- For a list of Svbonny-approved accessories for your item, visit the following website:
<http://www.Svbonny.com>

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EU Importer: Germany Retevis Technology GmbH
Address: Uetzenacker 29,38176 wendeburg

SVBONY

Warranty Card

Product Model

Purchasing Date

Defect Reason

Dealer Name

Telephone

User's Name

User's Address

User's Email

SYBONY

Warranty Card

Remarks:

1. This guarantee card should be kept by the user, no replacement if lost.
2. Most new products carry a one-year manufacturer's warranty from the date of purchase.
3. The user can get warranty and after-sales service as below:
 - Contact the seller where you buy.
4. For warranty service, you will need to provide a receipt proof of purchase from the actual seller for verification

Exclusions from Warranty Coverage:

1. To any product damaged by accident.
2. In the event of misuse or abuse of the product or as a result of unauthorized alterations or repairs.
3. If the serial number has been altered, defaced, or removed.